




### **Connecting Microscope Cameras HDCAM6B & HDCAM7 to a Windows Computer**


1. Download ToupView software.
2. Install software onto your computer. (Make sure the software is closed before connecting the camera.)
3. Connect the camera to your computer using the USB port, and turn on the camera.
4. Open the ToupView software, and the software should automatically detect your camera. If the software does not automatically detect your camera, follow the steps described in the subsequent pages to allow your computer to recognize the camera.



# How to use winusb driver on win8 and above system

1、 Find unrecognized cameras in Device Manager


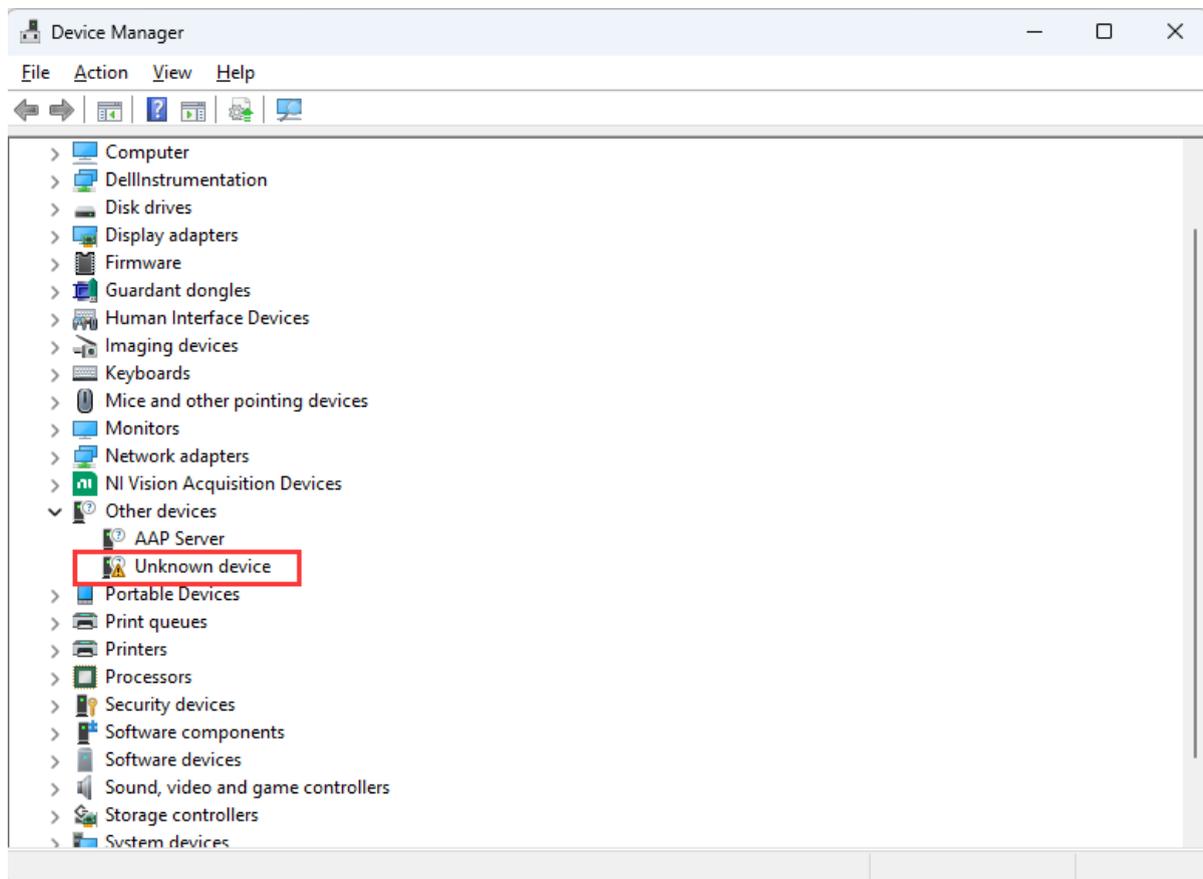


2、 Right-click to update the driver

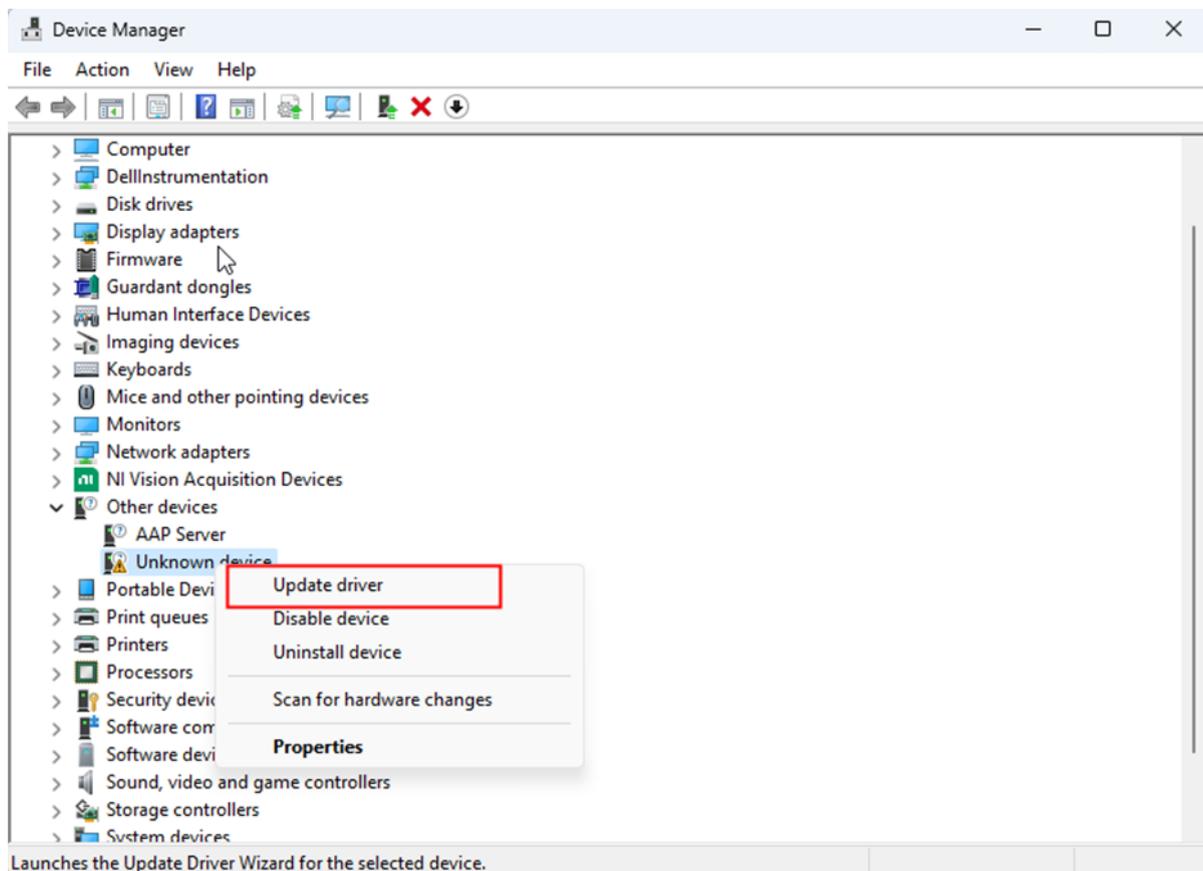


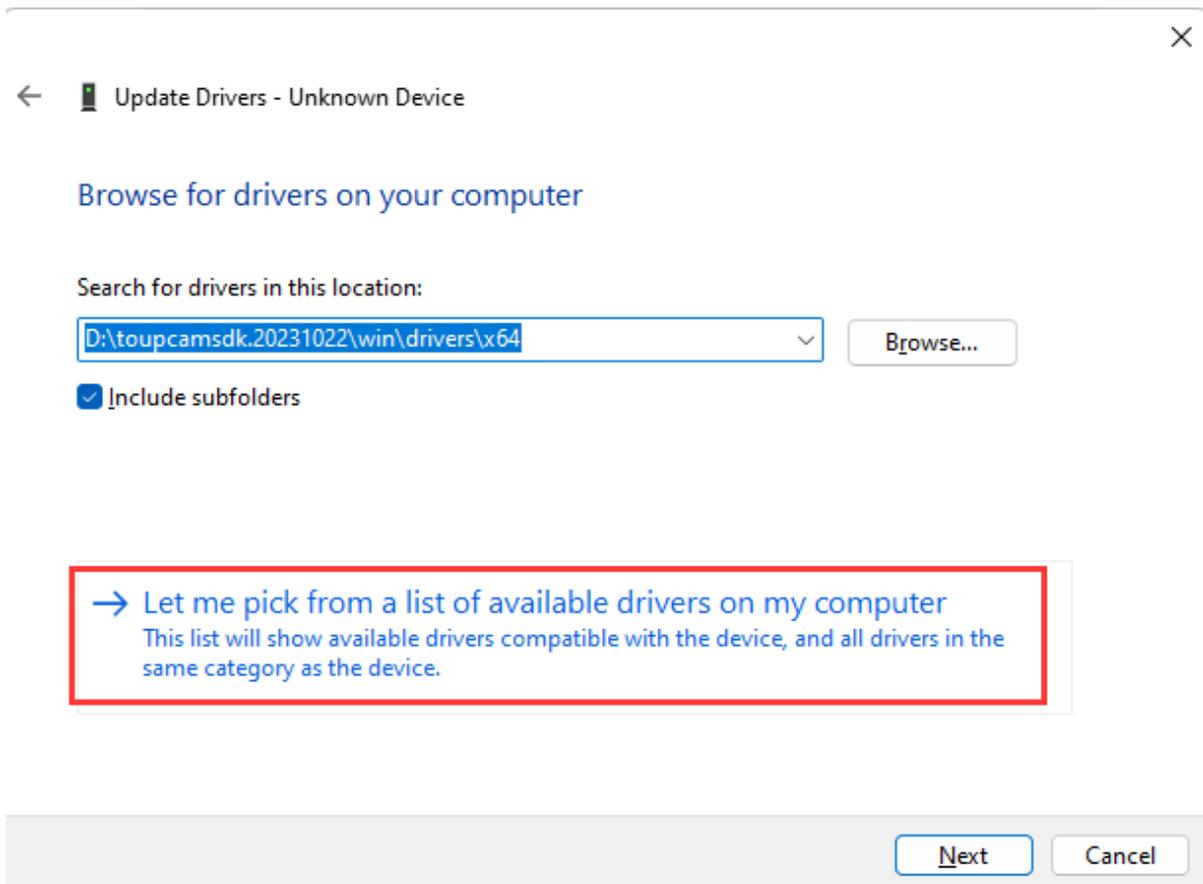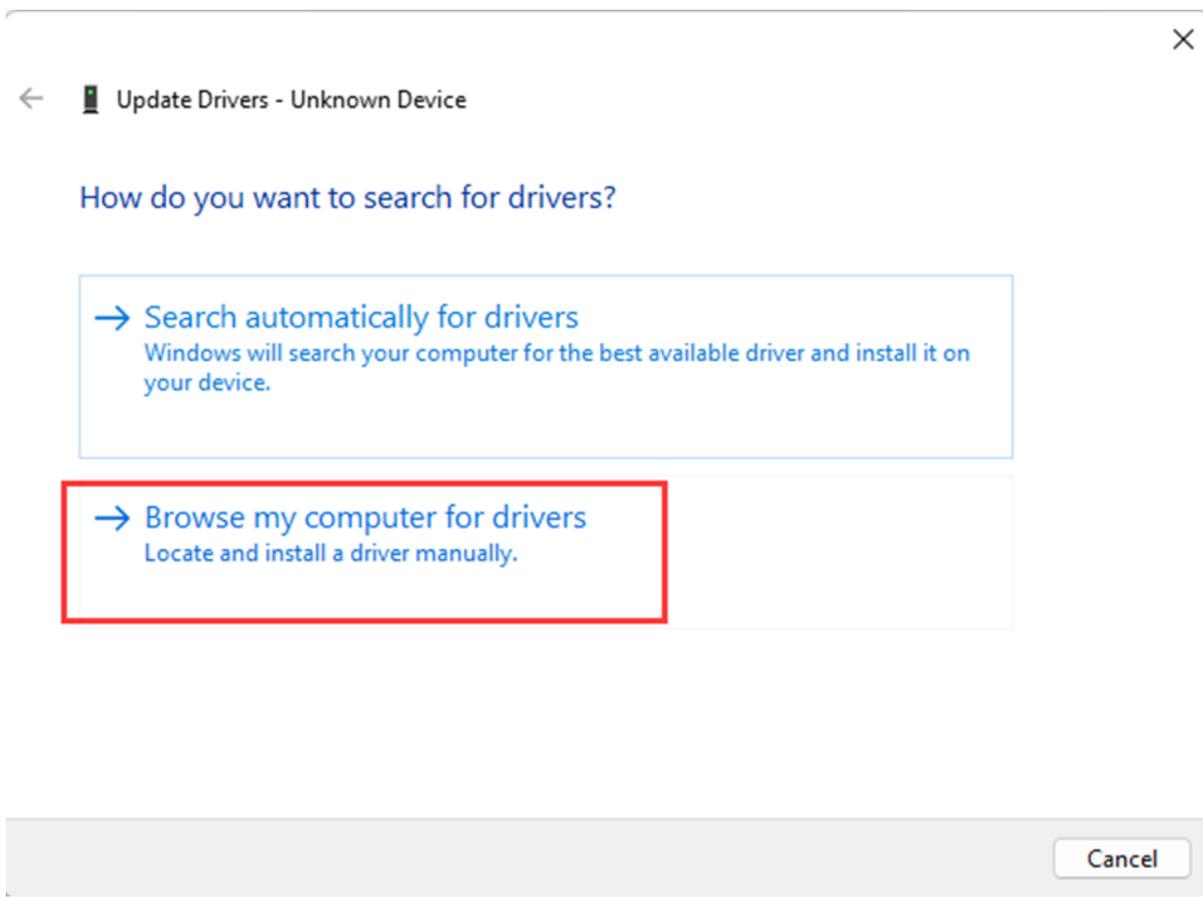

3、 Select the universal Serial Bus device

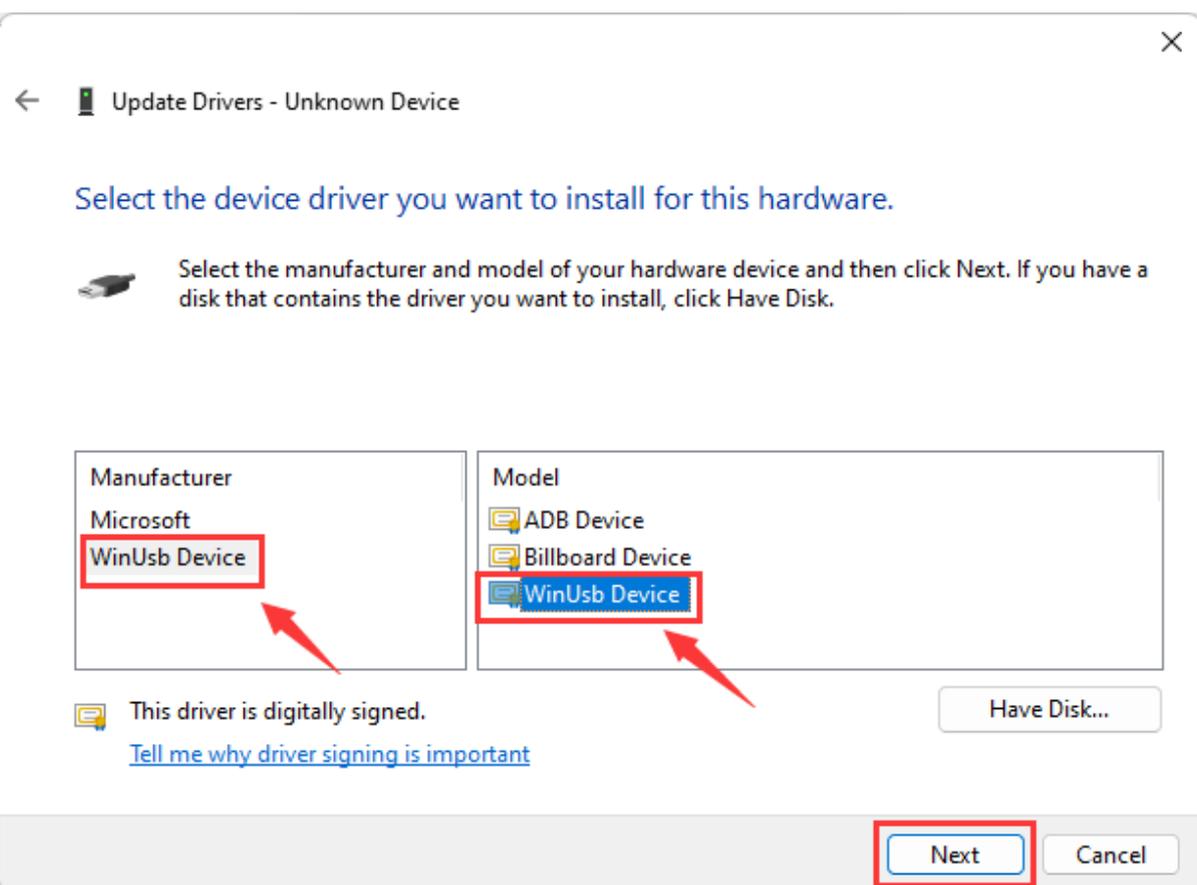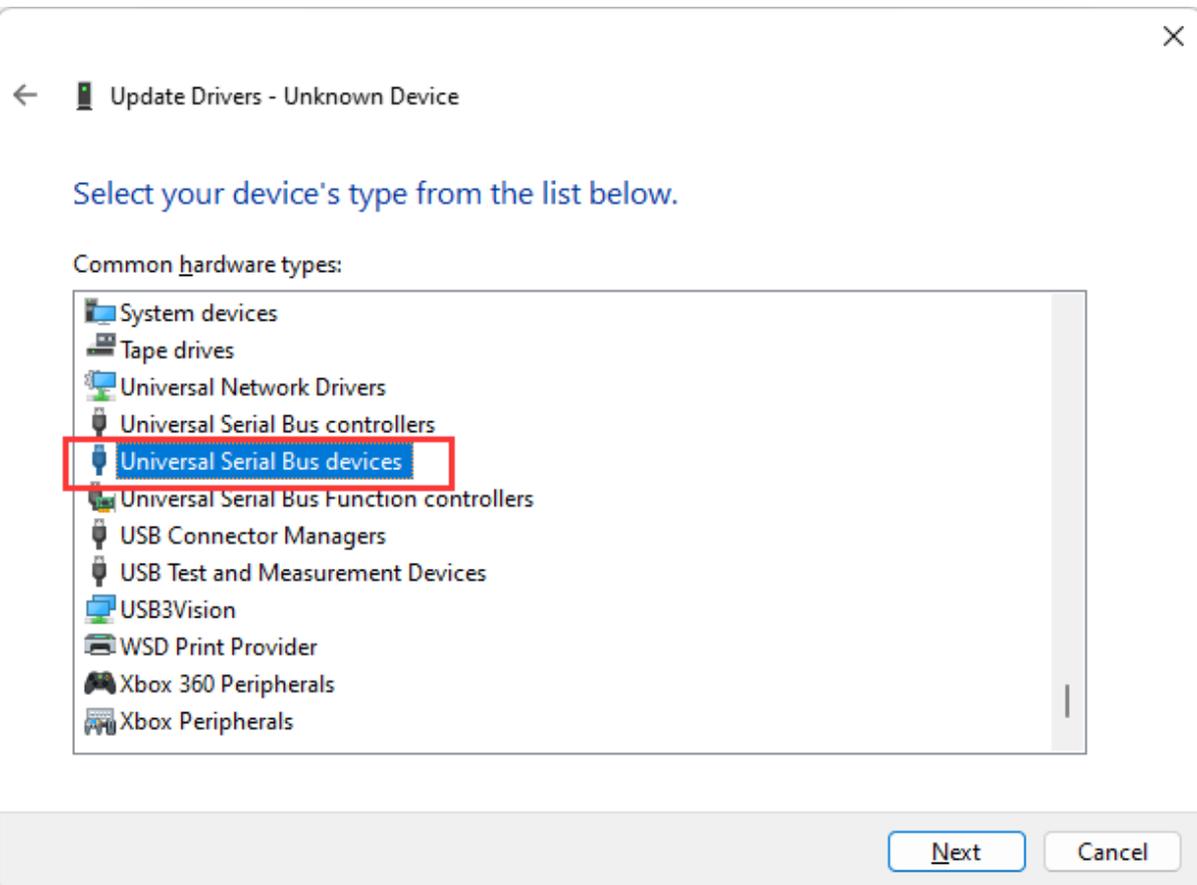

4. When a warning is displayed, click Yes to forcibly use the winusb driver

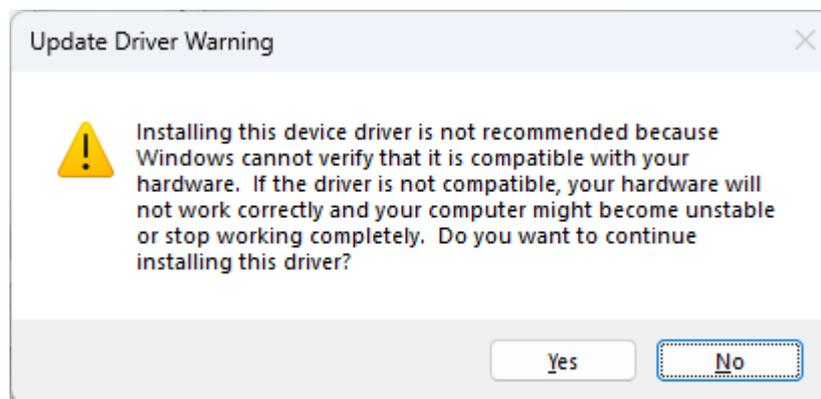

5. Finally, the camera is recognized as a winusb device in the Device Manager

